STERI
LIFT ®

CO,RA®

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The STERI LIFT is a drum, bin or bag positioner/lift/mixer that can be used in sterile rooms, and which allows quick and efficient handling.

The experience in facing the complexity of the strict norms in the pharmaceutical field and the ability of manufacturing stainless steel products has allowed us to introduce equipment suitable for the transportation and handling of products for chemical, pharmaceutical, cosmetic and food industries, marked by the continuous innovation of production processes into the market.

The main characteristics of the CO.RA. STERI LIFT are the hydraulic piston, completely enclosed in a casing. Any kind of contamination is avoided thanks to the perfect seal, as all movements are motorized through a motor/hydraulic pump with electric drive.

The motor is located inside the base of the STERI LIFT itself, while its casing (in AISI 304 stainless steel) provides insulation from the working environment.

The absence of exposed chains, gears, guides, eyelets and electrical and mechanical parts, as well as its realization with rounded surfaces, makes cleaning operations of the STERI LIFT easy.

The STERI LIFT complies with the recommendations of GMP (Good Manufacturing Practice) ensuring maximum protection of health and the prevention of accidents to the operator during the work cycle.

The surface of the not in contact stainless steel parts is satin finished Ra 0,8. Upon request and depending on the zone where it is installed, the Steri Lift can be supplied completely polished.

The hydraulic fluid consists of oil, or upon request, of hydraulic food grade liquid. Particular systems are foreseen to guarantee the perfect seal of the hydraulic fluid preventing it from getting in contact with the working environment in case of leaks or damage.

The STERI LIFT, once correctly installed in its service position, is always ready to be used. Its typical installation is on the ground, but special ceiling mounted versions are also available.

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Advantages:

- Simple and compact design
- Designed and manufactured based on the enduser's specifications.
- Easy to inspect and clean.
- Material: stainless steel with a silicone seal (certificates upon request).
- Manufactured in compliance with G.M.P.
- Easy and quick to maintenance.
- Adaptable to different kinds of containers.
- All system mechanisms completely encased.

Applications:

- Product handling to fill process machines.
- Container or Bin blending.

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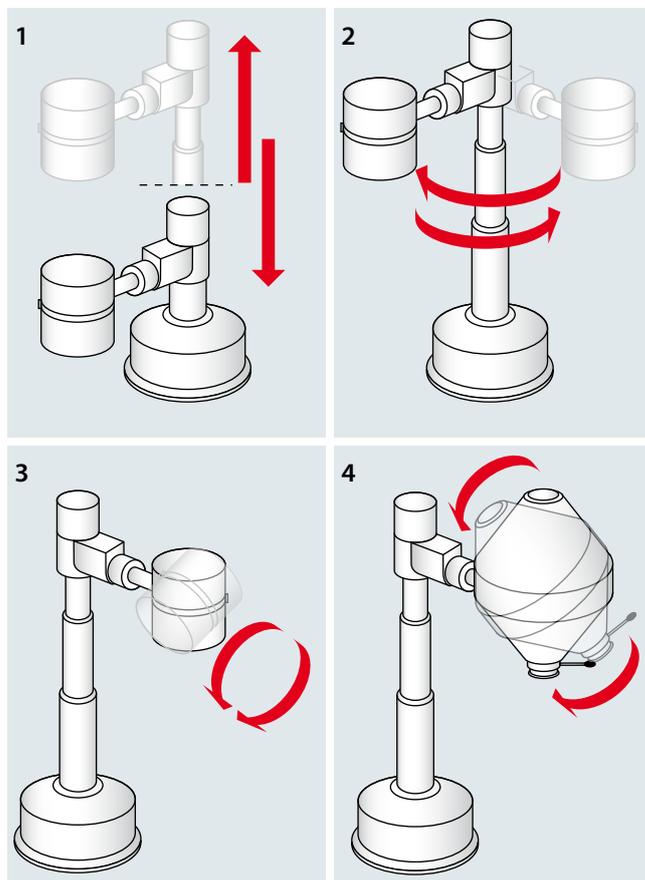
The STERI LIFT can perform three operations: manual (with the exception of the lifting/lowering movements), motorized (controlled by a pushbutton) or automatic (controlled by a PLC).

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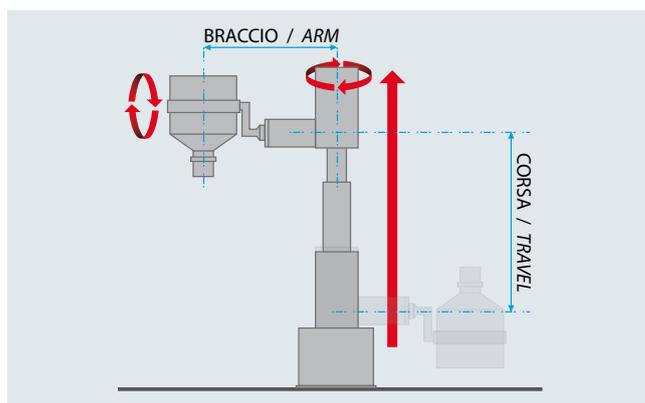
Based on the drive chosen by the customer the machine can be supplied with an automatic or manual work cycle. For the manual work cycle all the movements are controlled by the operator with a control keyboard, while for the automatic work cycle all the machine movements can be carried out with HMI panel or with the main switchboard. If a STERI LIFT is sold with an automatic work cycle, the machine is sold equipped with sensors or travel stops that limit the movements automatically.

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Attention: The type of drive for the various movements of the STERI LIFT is also chosen based on the type of drum/bin and on the mass to be handled by the machine.

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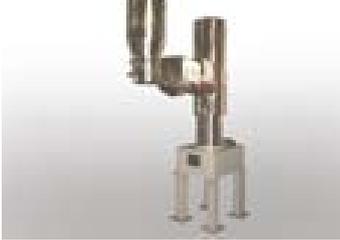
Base

Its function is to ensure the steadiness of the STERI LIFT as it is fixed to the ground or to the ceiling in the working environment. The top consists of a cover fixed with screws and can be lifted for maintenance operations.

All the hydraulic components that allow the movements of the machine, can be installed inside the base.

Below are listed the main types of STERI LIFT bases that can be installed depending on the working environment and the intended use of the machine.

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
Hood

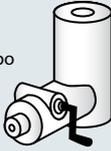The STERI LIFT hood is the structure fixed to the lifting piston. This element, which is joined to the lifting piston, transmits the movements of the latter to the Bin allowing the translation, capsiding, lifting, lowering and mixing movements.

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




Gripping System

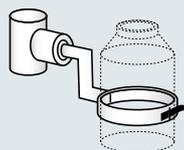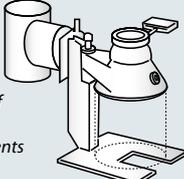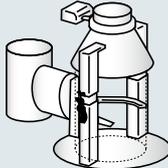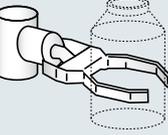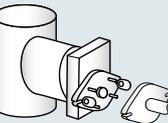The Gripping system is the element that locks the bin/drum to the STERI LIFT so it can be moved. The Gripping System is an accessory which is customized case by case according to the customer's needs.

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STERI LIFT + BICONO
STERI LIFT + BICONE

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Control Panel

All the buttons necessary for the operator to carry out all the movements of the machine are installed on the control panel of STERI LIFT. This kind of panel is installed in a such a position that prevents the operator from entering the working range of the machine during its movements. For security reason the control panel has been designed in such a way that it must be used with both hands.

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Electric Control Panel

The STERI LIFT control panels are designed and manufactured as an integral part of the machine and the complete knowledge of the manufacturing process allows us to define all the movements and the controls necessary for the operator to use the machine correctly and safely. All the buttons and lights that allow the operator to verify/check the status of the machine during all phases of the working cycle are positioned on the STERI LIFT control panel. In the operation of complex systems, where the STERI LIFT must be electrically interfaced with other system machines, the technology adopted by CO.RA. is PLC with touch screen for a simple and direct user interface. The use of the touch screen panels offers many advantages over traditional systems such as: system customization, extension of the functionality of the machine, setting of the working cycle, batch discharge batch handling, use of machine with a password and all the features required by the enduser.

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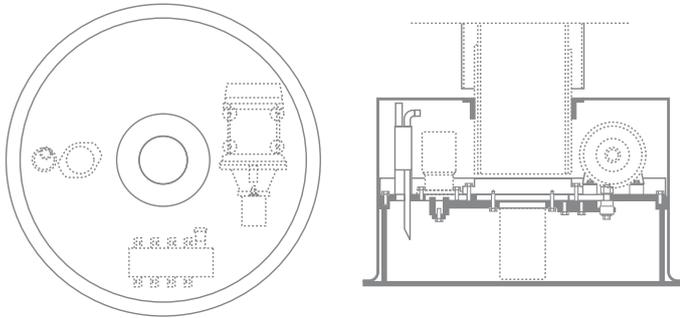
Hydraulic Unit

Tank where power equipment, control equipment of the lifting piston (motor/pump, control valves) are assembled. The hydraulic unit can be installed inside the base or, if it is possible, it can be separated and installed in a lining. This second option allows to reduce the dimensions of the base leaving more space for the movement of materials and operator inside the working environment.

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
Centralina idraulica a bordo
On board hydraulic unit

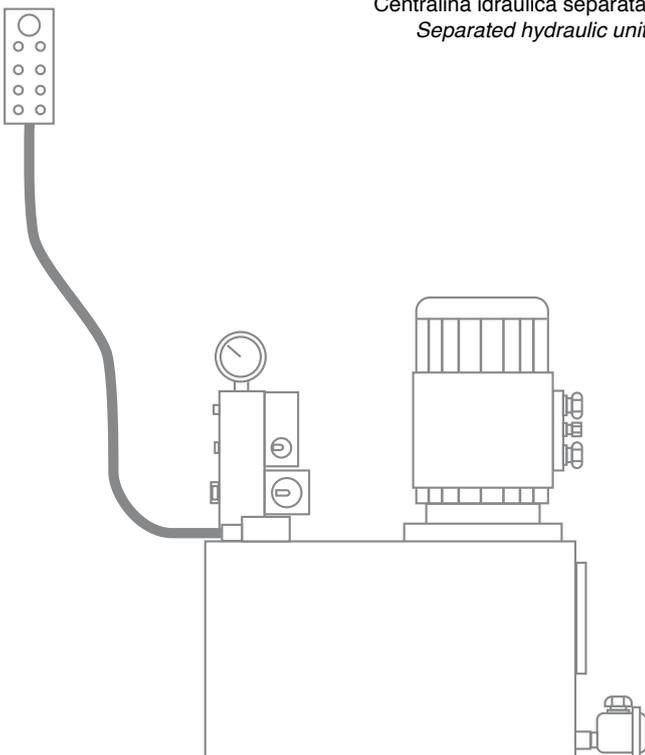
Centralina idraulica separata
Separated hydraulic unit

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
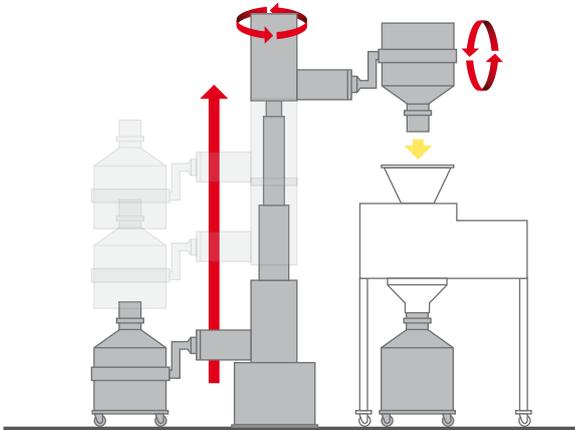
EXAMPLES OF STERI LIFT APPLICATIONS


Carico mulino - Mill feeding

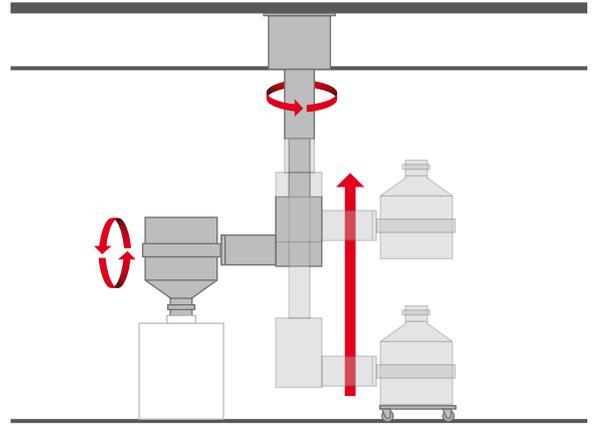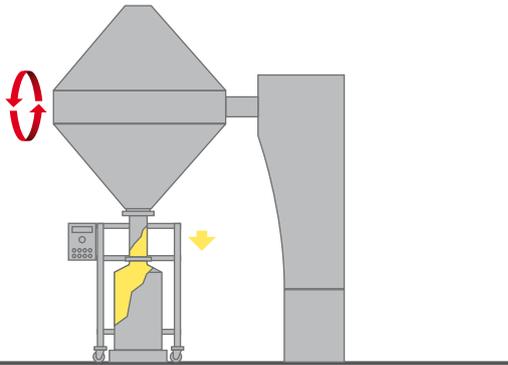
STERI LIFT appeso al soffitto - Ceiling mounted STERI LIFT

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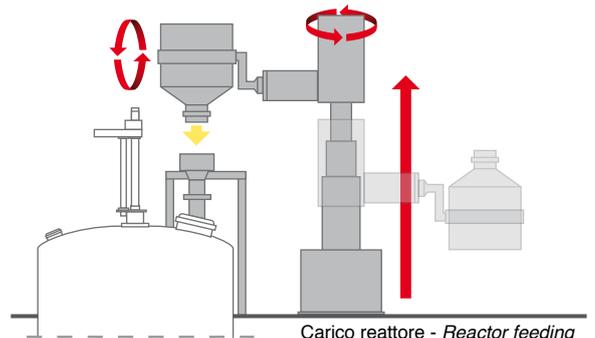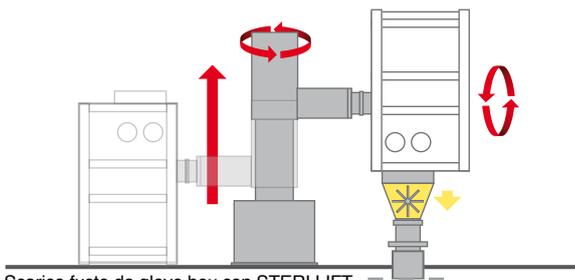
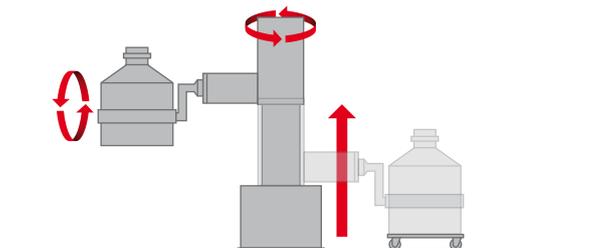
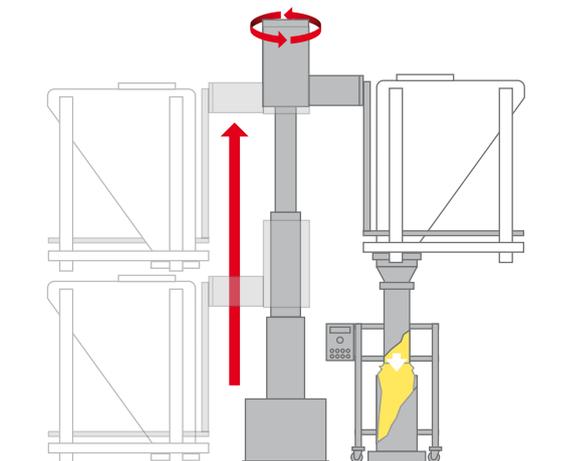
Miscelazione fusto con STERI LIFT - Drum mixing with STERI LIFT

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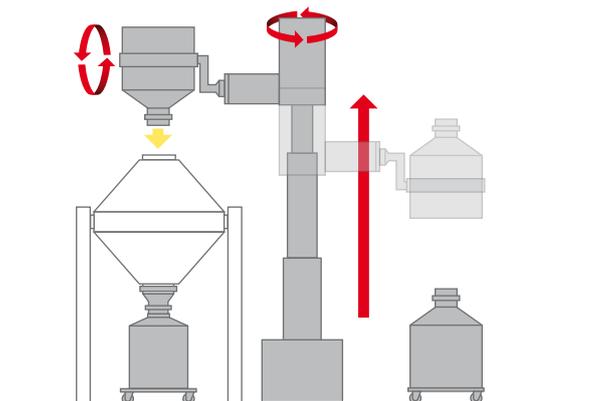SAFETY OF THE MACHINE

The STERI LIFT is designed to provide intrinsic safety to its structure, to prevent accidents or dangerous situations while using the machine.

- 1 In the event of a collision during the translation movement, the hood-gripping element group does not provoke any damage because the piston, which it is fixed to, consist of telescopically connected stems.*
- 2 The hydraulic unit is designed to maintain the STERI LIFT in the same position in case of malfunction thus avoiding crushing or collision with things or persons under or near the machine in that moment.*
- 3 To avoid that the operator may be accidentally bumped by the machine it is supplied with automatic working or mixing cycle, zone demarcation signals such as photocells, door lock sensors or laser scanners can be supplied. Once they are turned on they interrupt all the movements of the machine putting it in emergency mode.*

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OPTIONAL BENEFITS OF STERI LIFT

- *Adjustable product mixing speed*
- *Adjustable speed for an accurate match/connection to the inlet of the machine to serve*
- *Control valve to optimize the ascending and descending speed of the STERI LIFT*
- *Control the opening and closing of the product discharge valve by means of a keyboard.*

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ACCESSORIES

Besides designing and manufacturing the STERI LIFT CO.RA. designs a wide range of spare parts functional to the machine, such as:

- *Bins from 500 to 3000 Lt*
- *Containers*
- *Bi-Cones*
- *Hoppers*
- *Intercepting and dosing valves*
- *Roller conveyors for the handling of the container inside the STERI LIFT loading area*
- *Trolley for drum handling which are customizable to specific customer requirements.*

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EXAMPLE OF STERI LIFT ACCESSORIES

Bin con SOLE VALVE - Bin with SOLE VALVE

Bin 750 Lt con SOLE VALVE - 750 Lt Bin with SOLE VALVE

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DOCUMENTATION AND CERTIFICATES

The documentation supplied is:

- Material certificates of the parts in contact with the product
- O & M Manual
- Supply layout
- Electric and pneumatic diagrams
- P & ID
- CE certificate of conformity

Documentation which can be supplied on request:

- Functional specifications
- FAT e SAT protocols
- IQ ed OQ protocols
- Surface finish certificate for the parts in contact with the product

The STERI LIFT complies with the Machinery Directive, and also can be supplied for use in ATEX classified environments. The ATEX zone marking for the STERI LIFT is following:

II 2 G c k b 88°C (T5) 0°C Ta+40°C X
II 2 D c k b 70°C 0°C Ta+40°C X

Services related to the supply

The services that can be related to the sale of the machine are:

- Installation
- Start-up assistance of our specialized technicians
- Staff training
- Plan maintenance carried out by our technicians
- Assistance within 48 hours from your claim

The STERI LIFT is designed and manufactured respecting customer requirements, and to better guarantee this service it is possible to carry out tests in our facilities with one of the STERI LIFT is installed in our Technology Center.

USEFUL INFORMATION

For further information please contact our Sales Dept by phone +39 0583 20590 or by e-mail: sales@coraitaly.net

REG. No. 5762/2

REG. No. IT-71044

Atex 2014/34/UE

CO.RA. S.r.l. - 55011 Spianate - Altopascio (Lucca) - Italy - Loc. Chiappini, 51
Tel. +39 0583 20590 r.a. - Fax +39 0583 20481 - info@coraitaly.net - www.coraitaly.net