TABLET
VALVE

CO,RA[®]

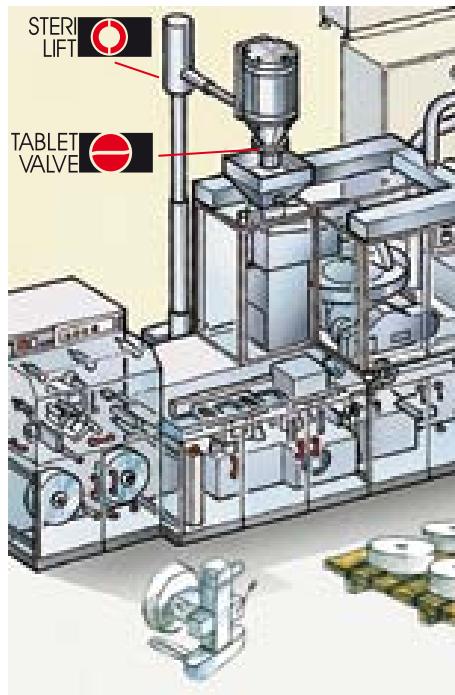
TABLET VALVE

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CO.RA. TABLET VALVE solves the problem of interception and monitoring of flows in a pharmaceutical plant and particularly allows handling of capsules, pills, tablets, rubber and silicone caps, using the same principle as the highly successful SOLE VALVE.

Its implementation derived from careful observation by CO.RA. of the operator's requirements during the most delicate pharmaceutical processes in a sterile room, especially the cleaning phase. The ergonomics of the parts of which it is composed make it easier for the operator during the disassembly and re-assembly of the valve.

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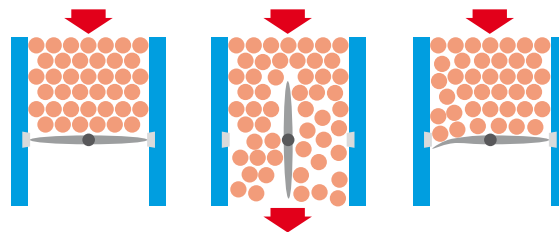
EQUIPMENT SERVED

- Packaging machines feeding
- Bottom valve for tablets, capsules, silicone caps, rubber caps containers
- More than 5000 different applications executed

The TABLET VALVE vane is coated and its outer surface is flexible. This allows opening/closing of the valve without damaging the product. The silicone is in accordance with FDA 21 CFR 177.2600 Standards.

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Product is held without any damage and can continue its flow as required.

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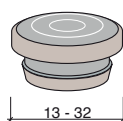
The speed and the stop motion are adjustable according to the requirements thanks to a control panel interfaced to the valve.


The range of half-bodies available and interchangeable, allows the user to make the application more suitable according to his disassembly needs.

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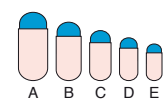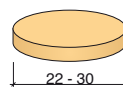
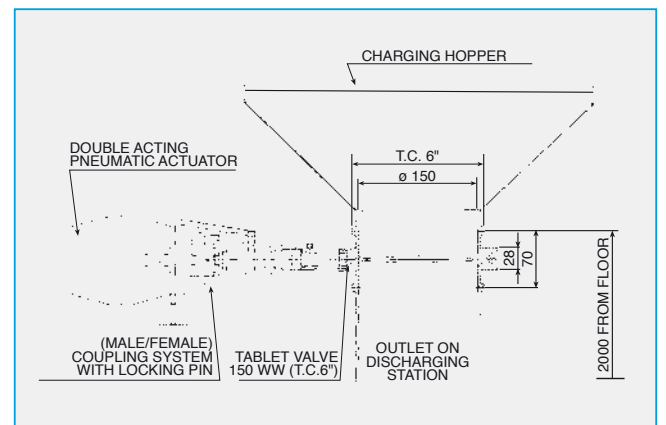
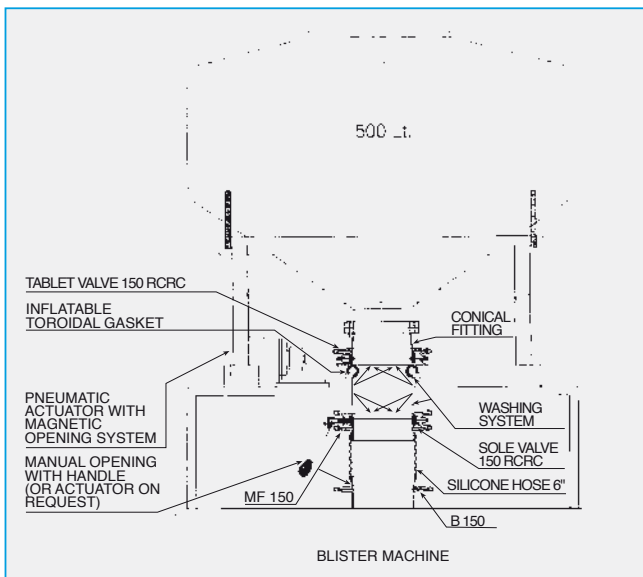
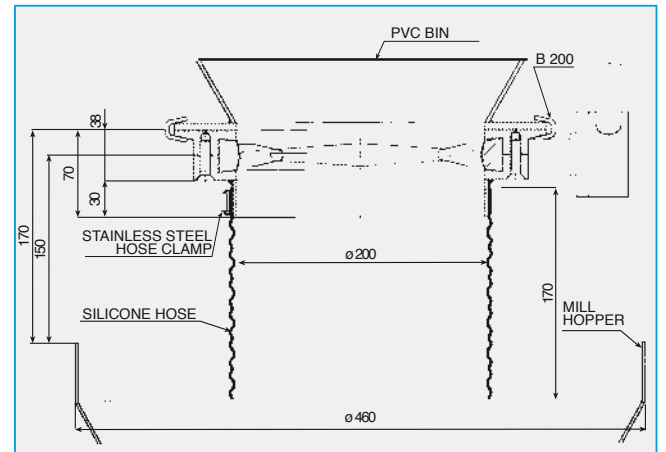
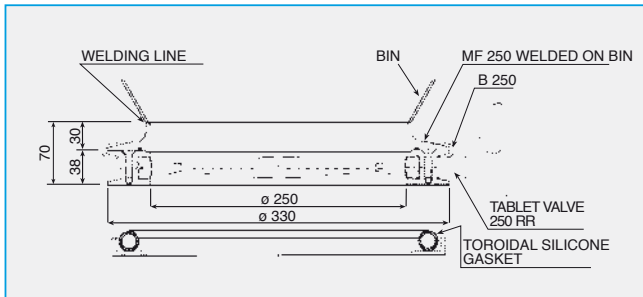
APPLICATIONS

The unique design of the TABLET VALVE provides a flow control system in less space and with smaller f.t.f. compared to conventional valves.

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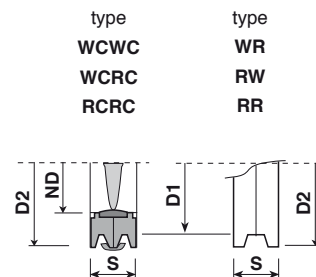
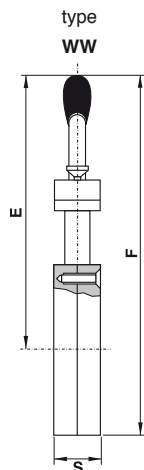
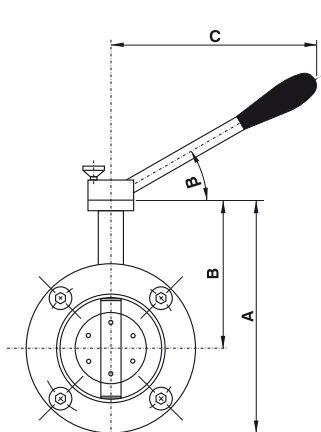
TECHNICAL FEATURES

- AISI 316L (DIN1.4404) satin finished body (mirror polished upon request).
- Silicone butterfly certified F.D.A. 21 CFR 177.2600.
- No O-Ring (OR) on the shaft.
- Easy to disassemble, inspect, clean, particularly the "quick release" series with "CLAMP".
- Reduced weight and dimensions.
- Materials: stainless steel and silicone (certificates upon request).
- Manufactured according to G.M.P.
- Suitable for tablets, pills and granules.
- Traceability over time (batch number).
- Interceptor element interchangeable with ROTARY VALVE e SOLE VALVE.
- Each valve bears the CO.RA. trademark and identification batch number (ex: CO.RA. I/00/298/GP).
- Anodized aluminium pneumatic actuator.
- Air command 2 ÷ 6 bar (30 ÷ 87 psi).
- Consumption per drive, pression 5.6 bar (80 psi): 3AD=1.68 Lt/m.

Advantages:

- Size and weight reduced, compared to traditional valves.
- No lubricants.
- Sterilizable also in autoclave.

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


Stainless steel handle on/off, different positions with bachelite knob (silicone knob upon request)

- Upon request it can be supplied:
- 45° handle
 - Quick release Sanitary handle
 - Removable security handle

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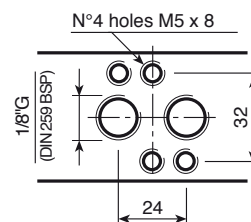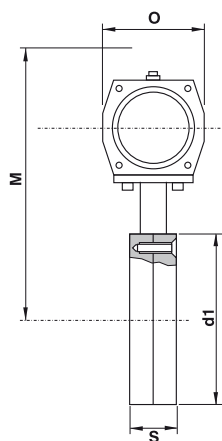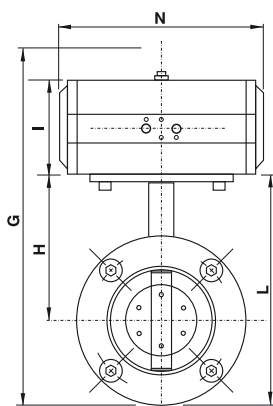*Double acting actuator (AD) anodized aluminum to ISO 5211
 Air supply: 4 ÷ 6 bar (58 ÷ 87 psi)
 Air consumption to 5,6 bar (80 psi)
 AD 63-180 = about 0.26 liters
 AD 85-180 = about 0.55 liters*

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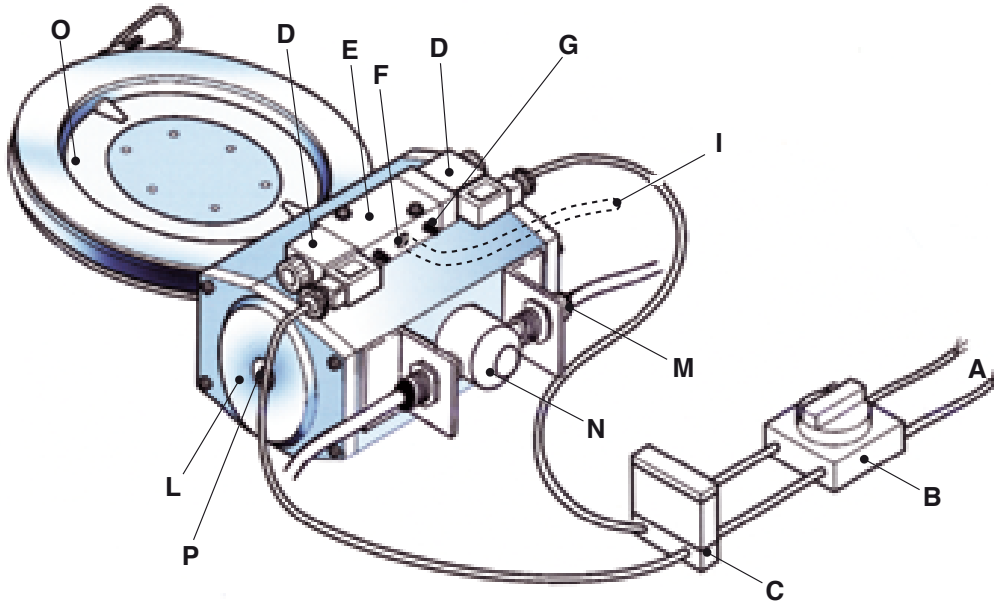
The mounting flange can be supplied with male end MF/M if connected to valves and female end MF/F with OR and MF/F with OR; the same for blanking plate P, P/M, P/F, P/F are made of silicone if requested of other elastomers.

*Ask for the list of all hand levers, supports and opening systems.

**Also other types of coupling available: Bayonet, chute, etc.

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
Example of assembly plan with double acting (AD) pneumatic actuator

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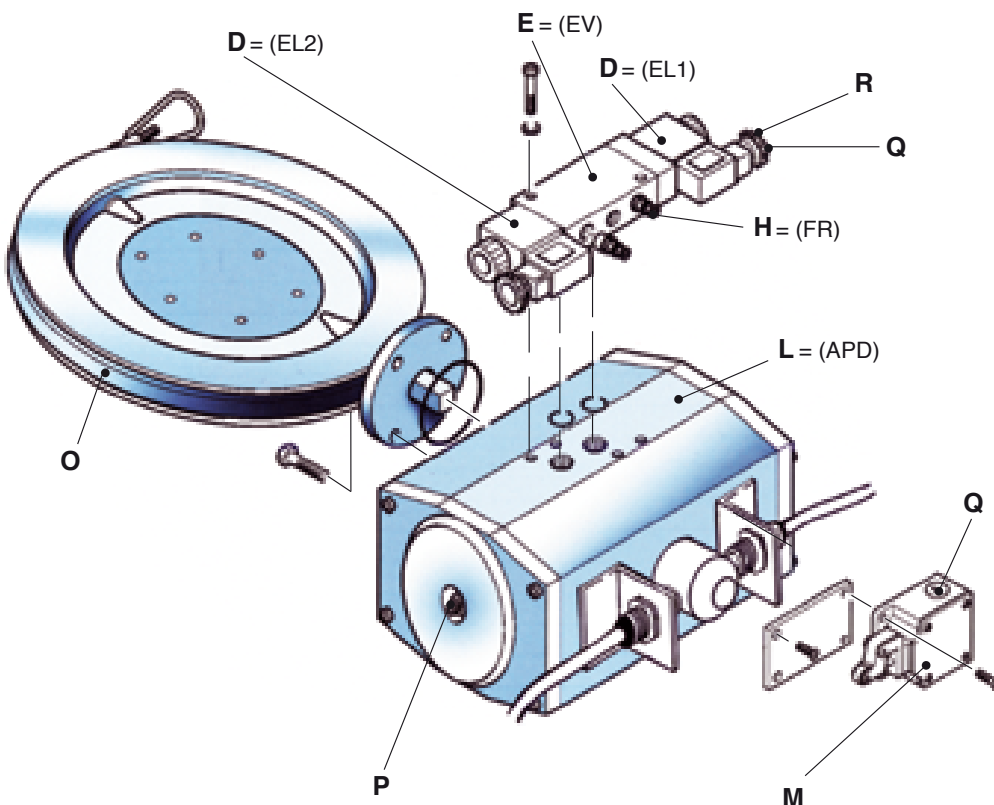
Exploded view for spare parts ordering (for spares for each main component ask for detail drawings)

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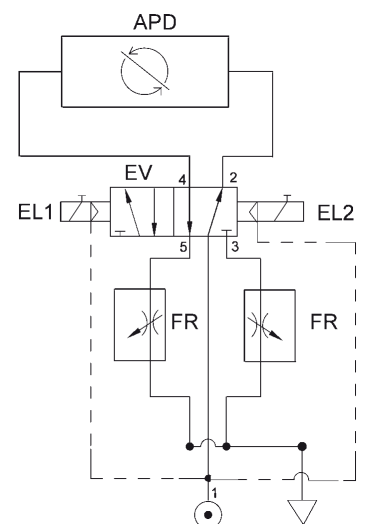**TABLET VALVE heavy "H"
series stainless steel – silicone**

Body made from solid

Suitable for:

- Flow control of very hard pills, granules
- Flow regulation of hard, small tablets

The stainless steel halfbodies have no gasket.

Actuation: by handle or pneumatic actuator.

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**TABLET VALVE "K"series
plastic – silicone
(version available for all types)**

It is designed to be used when the stainless steel body is not usable or required.

Useful to be mounted on PVC bin. Available without gasket upon request.

Actuation: handle or with pneumatic actuator.

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Comando: solo con leva.

**Economic TABLET VALVE "E"
serie, stainless steel – silicone
(available only model "U" type)**

Very light, obtained from AISI 316 pipe (DIN 1.4401), very economic.

Suitable for:

- flow shut-off of granules, empty capsules, tablets, pills.
- control of air conditioning in the bearings (cleanroom, air conditioned rooms, sterile room air conducts).
- regulation of very hot air in sterile room conducted drying ovens, drying tunnel.
- Quick release with special «CLAMP» or welded directly on the piping.

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DOCUMENTATION AND CERTIFICATES

The documentation that can be supplied with the valve is:

- Use and maintenance manual
- Material certificates of the parts in contact with product
- Certificate of surface roughness for the parts in contact and not in contact with product
- Functional specifications
- Certificate of conformity CE
- FAT and SAT protocols
- IQ and OQ protocols

The Tablet Valve complies with Machinery Directive, is CE marked and also can be supplied for Atex zones.

The valve markings can be as follows:

II 1GD/2GD c 85°C (T6) 0°C Ta+40°

II 1D/2GD c 85°C (T6) 0°C Ta+40°

The Tablet Valve to be marked / used in Atex environment 1GD/2GD must have the butterfly made of black conductive silicone (F.D.A. approved). For the ATEX marking 1D/2GD contact our sales department to know the available elastomers (all F.D.A. approved)

Services related to the supply:

The services that can be related to the sale of the supply are:

- Assistance within 48 hours from malfunctioning.

The TABLET VALVE is designed and manufactured respecting customer requirements, and to better guarantee this service it is possible to make test at our premises with one of our TABLET VALVE in our Technology Center.

USEFUL INFORMATION

We suggest you to use the above indicated abbreviations (W, WC, WR, WCRC, ...) to easy fill in your P.O.; please consider the halfbody are usually made of AISI 316L. The seal is elastomer.

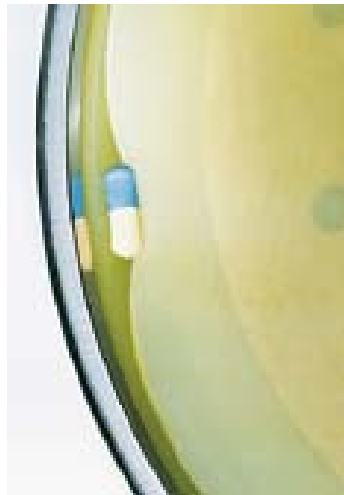
Each valve with same diameter has interchangeability about half bodies, supports, rotors, seals, actioning, ecc.

The clamps are the same for quick clean and quick release.

Gasket available in several material and types are in accordance with FDA, for more information please ask for SOLE VALVE catalogues.

For further information please contact our Sales Dept by phone +39 0583 20590 or by e-mail: sales@coraitaly.net

Reproduction of part all of the contents of this publication without the written consent of CO.RA. s.r.l. is forbidden

REG. No. 5762/2

REG. No. IT-71044

Atex 2014/34/UE

CO.RA. S.r.l. - 55011 Spianate - Altopascio (Lucca) - Italy - Loc. Chiappini, 51
Tel. +39 0583 20590 r.a. - Fax +39 0583 20481 - info@coraitaly.net - www.coraitaly.net